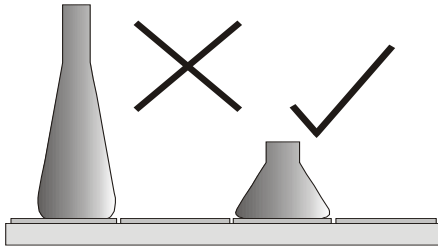
Always use protective goggles and gloves while handling glassware!

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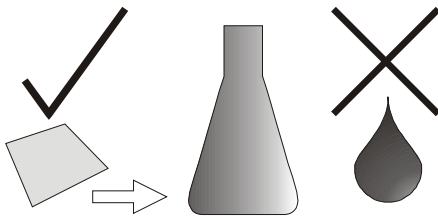
For containers with a wide, flat base only. Large Erlenmeyer flasks (e.g. 3000 mL) will stick stronger than small ones (e.g. 100 mL).

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Before putting on containers:

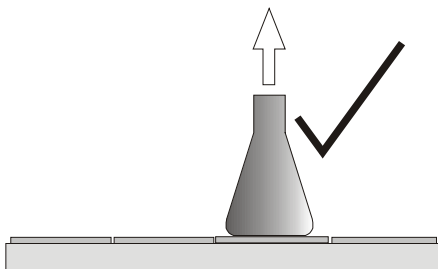
- Check that bottom and adhesive matting is dry and clean
- Check flasks for any damages, **never use damaged flasks!**

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Tug gently to ensure each flask is stuck. Keep in mind that formation of condense water might occur at low temperatures or using the timer function, possibly causing accidental release of containers.

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
To remove flask: Push or pull neck gently and evenly and wait a few seconds! With large containers, it can take up to 30 seconds until flask is released. Do not use much power!

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
Directions for use | Adhesive Matting „Sticky Stuff“

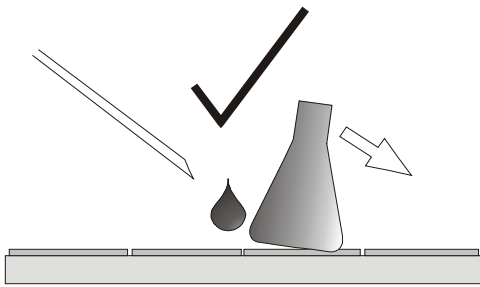Flasks which are firmly attached can be released with water. For containers with large bottom surface the adhesive matting may be covered partly with the protective foil supplied with each new tray.

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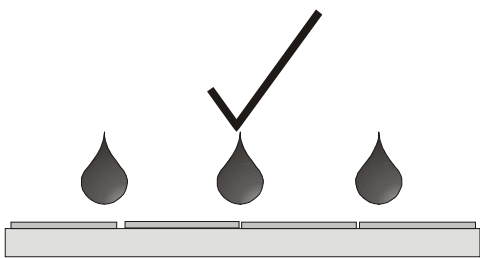
Clean regularly with water and mild detergent. Shrub vigorously with a brush or hard sponge and let dry over night. Do NOT use solvents! Disinfection with 70% ethanol or usual disinfectants, rinse thoroughly with water. If disinfection is done routinely, it might be necessary to replace the Sticky Stuff earlier.

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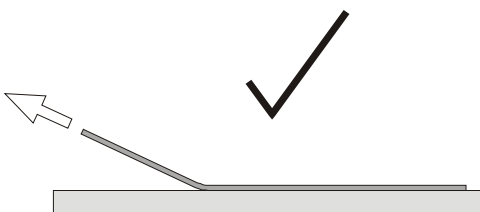
Changing of Adhesive Matting: Release Adhesive Matting on one side and pull obliquely away from the tray. Degrease tray with Acetone. Apply new patches with water (see installation instructions). Do not remove protective foil until first use.

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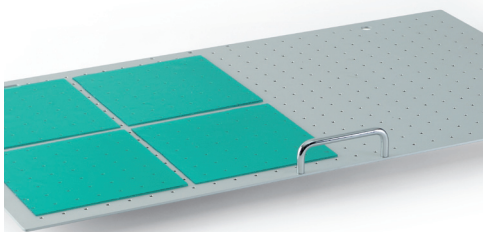It is possible to apply the Green Adhesive Matting to universal trays.

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Maximum shaking speeds:

Only valid for original Schott Duran® Erlenmeyer flasks with 20% filling volume and paper- or cotton stoppers. Glass and Adhesive matting must be totally undamaged, clean, dry and free of grease. All information supplied without liability.

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